



QUICK DECON COMPLETE POWDER MIX

ACTINIDE + TRANSITION METAL

DO NOT MIX WITH ACIDS /MAY RELEASE NOXIOUS GAS

Directions:

1. Completely dissolve powder in an appropriate container using clean (DI preferred) water.
2. Stir or shake the container to ensure adequate mixing
3. Apply a generous amount of solution to the surface to be decontaminated.
4. Let stand for up to 5 minutes.
5. Remove solution and contaminants with toweling or rinse and dispose of as radioactive waste.
6. Follow with cold water rinse and wipe to remove any residual solution.
7. The rinse effluent and/or wipes must be treated as radioactive waste.



Generally, not hazardous in normal handling;
however, good laboratory practices should be used.
Avoid long term exposure to skin or by inhalation.



www.rtlasersafety.com

770-332-0092

contact@rtlasersafety.com